



AtOMS Software Suite

User Manual

Release 1.2.1

Copyright AccuStrata 2021



Contents

Safety Symbols and Terms	2
Ultraviolet light warning.....	2
High voltage warning	2
Terms	2
Hardware Setup	3
Back Panel.....	3
Front Panel	3
HCL Compartment.....	3
External Hardware	4
Getting Started	4
Main Panel	6
Lamp Configuration Panel	7
File Saving Structure	7
Spectrum Tab.....	8
Peak Tracking Tab.....	9
Peak Selection Tab	10
Data Viewer Tab	11
Calibration Tab.....	12
System Menu	13
Setup Spectrometer Display	13
Sensors to Display.....	13
Edit Peak Description	13
List of Currently Defined Peaks	13
Peak Definition Creator	14
Hollow Cathode Settings	14
Lamp Current	15
Detector Menu.....	16
Set Integration Time	16
Set Scans to Average	16
Set Shutter Ratio	16
Select Chambers	17
Serial Communication	17
Hardware Setup.....	17
Software Setup	18
Communication Protocols.....	18
Outputs.....	18
Inputs.....	18
About.....	19
Contact.....	19

Safety Symbols and Terms

Ultraviolet light warning

Hollow cathode lamp light sources produce ultraviolet radiation. It is important to wear proper eye protection when adjusting the focusing optics within the system. The system should be powered down when exchanging hollow cathode lamps and the access door should be closed during operation.



High voltage warning

Hollow cathode lamps use high voltage power supplies which can pose an electrical hazard. Do not open the AtOMS box or break the tamper resistant seals located on the paneling. Keep the system away from other high-power electronics.



Terms

! Recommendation

Recommendations provide users information that may not apply to all use-cases but can generally be useful to know.

i Note

Notes are important information that should be kept in mind while proceeding.


! WARNING

Warnings are important safety information that must be followed during the operation of the AtOMS system.



Hardware Setup


Back Panel

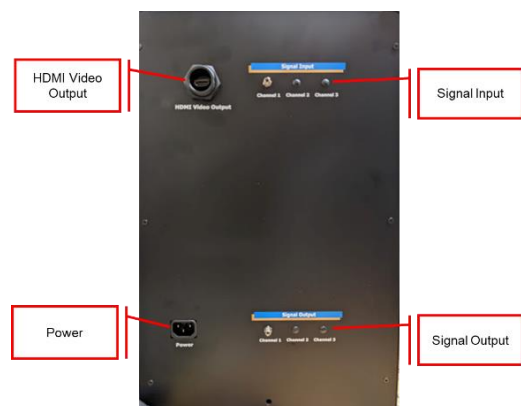

1. **Signal Output:** Light is emitted through these channels from the HCL. An SMA fiber optic should be attached leading to the emission optics on the system being monitored.
2. **Signal Input:** Return light coming through the stem being monitored comes into this SMA connector. An SMA fiber optic should be attached leading to the receiving optics on the system being monitored.
3. **HDMI Video Output:** Standard size HDMI video output for attaching to a display.
4. **Power:** AC Power Input 100-120V/5A

Front Panel

1. **USB 1+2:** USB A ports for attaching a keyboard and mouse or a USB drive.
2. **UV Light Indicator:** When the HCL is in use this light will be illuminated indicating the presence of UV light. Ensure proper eye protection if the HCL compartment needs to be opened while the indicator light is on.
3. **Power Button:** When the system is on the button will be illuminated in blue. To turn the system on press the button down to begin startup. To turn system off either shut down from Windows or press and hold Power Button for several seconds until system initiates safe shutdown.
4. **HCL Compartment:** This compartment is where the HCL and associated optics are located. While system is in use this compartment should always be shut and locked to ensure that UV light is not being released into the room.

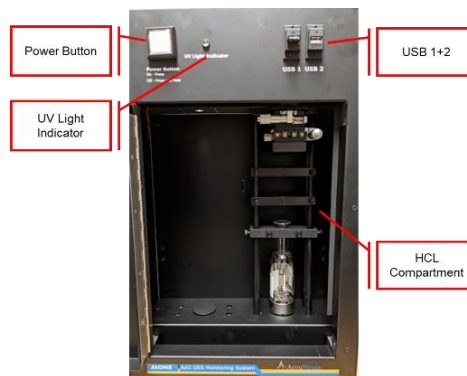

⚠ WARNING

This compartment is where the HCL and associated optics are located. While system is in use this compartment should always be shut and locked to ensure that UV light is not being released into the room.

HCL Compartment


1. **HCL Power Slot:** The system is designed for a standard 2-prong HCL. The HCL and connector are keyed so that the HCL is only capable of going in one direction.

⚠ WARNING

Ensure the AtOMS system is powered down when replacing Hollow Cathode Lamps.



2. **HCL Clip:** Clip to stabilize and hold the HCL in place. Clip is keyed in one direction and held in place by tightening the two hand screws.
3. **Slide Rail Optics:** Optics to focus the HCL onto the collector. Once maximum intensity has been reached tighten optics in place using the hand screws.
4. **XY Stage Adjustment:** Adjust XY stage in order to maximize light into the system when replacing HCL.
5. **Reference Power Adjustment:** Mechanical wheel used to adjust reference signal in software. It is recommended to run the system with the *initial reference and chamber signals within 5% of each other.*

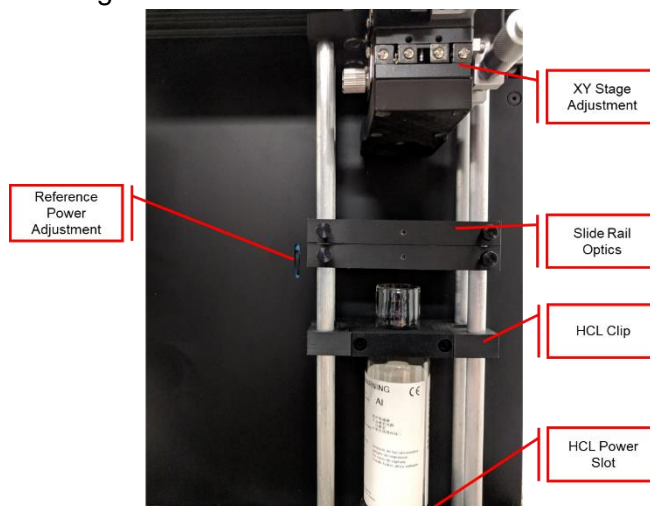

External Hardware

1. **Fiber Patch Cords:** For the standard system with 1 output and 1 input channel there will be 2 fiber patch cords to connect. The first patch cord will connect between the **Signal Output** on the AtOMS box and the **Emitter Optics** on the chamber. The second patch cord will connect between **Signal Input** on the AtOMS box and the **Collector Optics** on the chamber. To connect the patch cords, insert the SMA ferrule on each side into the SMA barrel connector and turn the threads clockwise until tight.
2. **Emitter Optics:** The optics used generally change based on a use-case basis. In most instances there will be an XY stage on the side of the emitter optics. This XY stage can be used to adjust the position of the optics in both directions using the spectrum tab on the software to obtain maximum signal throughput. The optics themselves are held in place via a 4-40 set screw.
3. **Collector Optics:** The collector optics will generally be attached to the chamber via a fixed position mount with a 4-40 set screw holding the optics to the mount. To remove the optics, loosen set screw before pulling collector optics out of the mount. In some cases, there will be an additional XY stage for the collector optics, in this case both the Emitter and Collector Optics XY stages will need to be positioned simultaneously to maximize signal throughput.

Getting Started

1. Plug system into 110V AC power source using provided power cord.
2. In HCL Compartment, plug desired HCL into HCL Power Slot and tighten in place using the HCL clip.
3. Turn on system using Power Button on front panel.
4. Confirm that the Fiber Patch Cords are connected to the Emitter and Collector Optics on the chamber
5. Once system has completed boot-up open AtOMS Software Suite



6. Set based software settings:

- a. Go to **System->Hollow Cathode Settings** and set the current provided to the HCL based on the specifications provided by the HCL manufacturer.

! WARNING


It is extremely important to stay within the current bounds that the manufacturer specifies as going outside of these bounds can damage the HCL.

- b. In software go to **Detector->Integration Settings** and set the value to 30ms.
 - c. In software go to **Detector->Scans to Average** and set the value to 5
 - d. In software go to **Detector->Set Shutter Ratio** and verify the value is set to 1.
7. Turn on HCL and begin taking data using Start Acquisition.
8. Optical Alignment
- a. Open the AccuAlign Software

! WARNING

AtOMS Software Suite must be **CLOSED** before starting AccuAlign Software.

- b. Select the chamber on the bottom of the window that will be optimized.
- c. Next, use the XY stages and goniometer on the chamber to maximize the signal through the chamber.
- d. Select the Internal button on the bottom of the window. In the HCL compartment use the Reference Power Adjustment wheel to make the primary peak for the HCL Reference and AAS+AES approximately equal. **You want these values to be within 5-10% of each other for best signal to noise.**

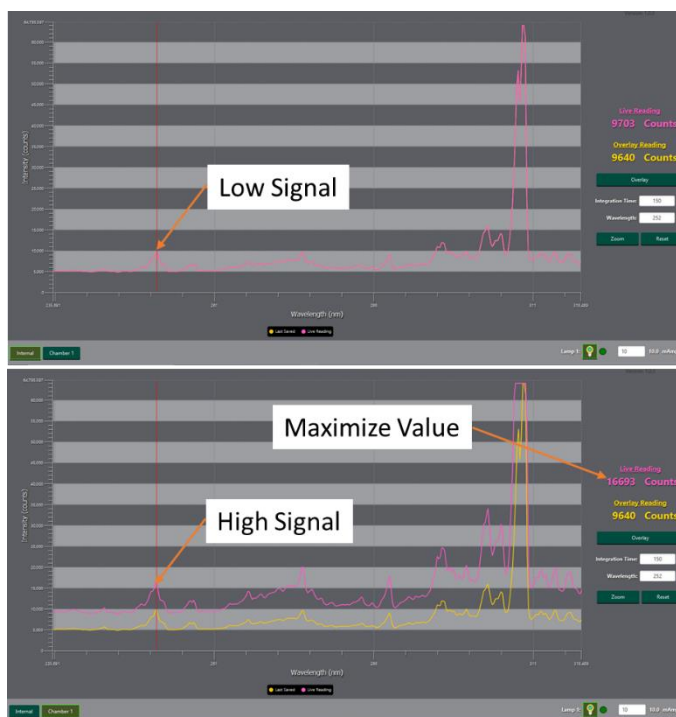

- e. Close the AccuAlign software and open the AtOMS Software Suite.

9. In **Detector->Integration Settings** change the integration time so that the AAS+AES and HCL Reference peak values are approximately 45,000 with neither value being greater than 50,000.



10. Once the integration time is determined adjust the *Scans to Average* to the desired level of signal to noise vs data collection time. The data collection time for a full scan in seconds is approximately equal to $4 \times (\text{Scans to Average}) \times (\text{Integration Time} / 1000)$
 11. Verify there is no deposition going on in the monitored chamber and take a reference.
-

Main Panel



Take Reference

To properly calculate absorption values, the AtOMS requires a reference reading. This reference should be taken when there is no deposition occurring. The reference should also be taken prior each run. Taking a reference will also automatically start acquiring data.



Start Data Acquisition

This button will start the data acquisition for the system.



Stop Data Acquisition

This button will stop the data acquisition for the system.



Start Recording

This button will start recording data that is being acquired in a new file. Recording can only occur if a reference has been taken.



Stop Recording

This button will stop recording data into the file that was created.



Lamp Control Display

This panel controls the hollow cathode lamp. The panel will auto-populate with the correct number of lamps that the system has installed. The circle indicates the current lamp status. The light bulb toggles the lamp.



Lamp Configuration Panel



Lamp Toggled On

Indicates that the user intends for the lamp to be on.



Lamp Toggled Off

Indicates that the user intends for the lamp to be off.



System Status-Lamp On

The AtOMS system is providing power to the Hollow Cathode Lamp.



System Status-Lamp Off


The AtOMS system is not providing power to the Hollow Cathode Lamp.

File Saving Structure

The AtOMS uses a Default file location of:

C:/Users/AtOMS/Documents/AAS Data

After the “Stop Recording” button  is pressed, a popup will open.

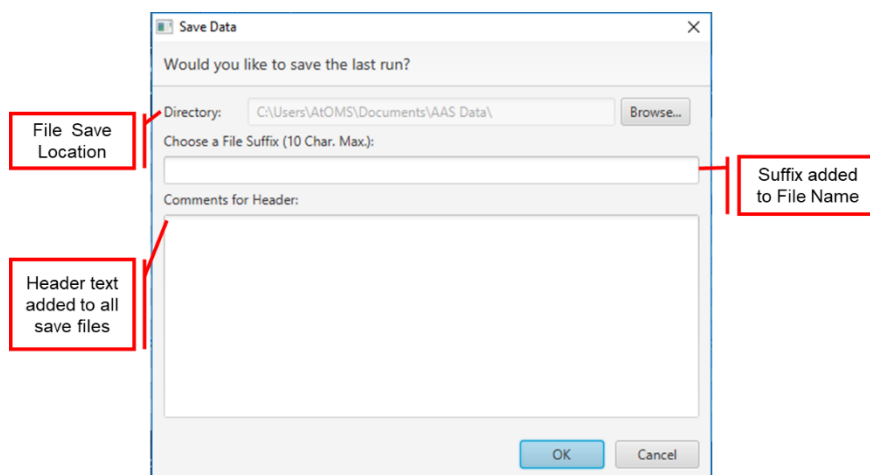

The **Directory** can be changed with the first “Directory” field. The **File Suffix** is text that will be appended to all files that are created. The **Header** field is appended inside every file that is saved.

The file Naming scheme is as follows:



AAS Data/YYYY_MM_DD/HH_MM_SS Suffix/

Within the folder there will be 4 .csv files and another folder that contains all processed data. The 4 .csv files are the raw data taken by the system. The processed data folder contains a file for each tracked element.

Spectrum Tab

The Spectrometer Tab Displays the raw spectral data that the AtOMS system is measuring. There are 4 types of data that can be displayed on this panel.

AAS + AES

Displays the total raw counts that have been measured. The signal will inherently contain both atomic absorption and atomic emission signal for each spectral line.

AES

Displays the atomic emission. This is taken by blocking the Hollow Cathode Lamp light and measuring the native emission from the chamber.

HCL Reference

Displays the “power measurement” of the Hollow Cathode Tube. This is taken by passing the light from the Hollow Cathode Tube directly to the spectrometer, bypassing the deposition chamber.

Baseline


Displays the baseline of the spectrometer. It is acquired by measuring the signal that is being received with the Hollow Cathode Lamp turned off and by passing the chamber.

Reference Measurement

Displays the total power of the Hollow Cathode Lamp that was taken during the reference phase of the acquisition.



Peak Tracking Tab

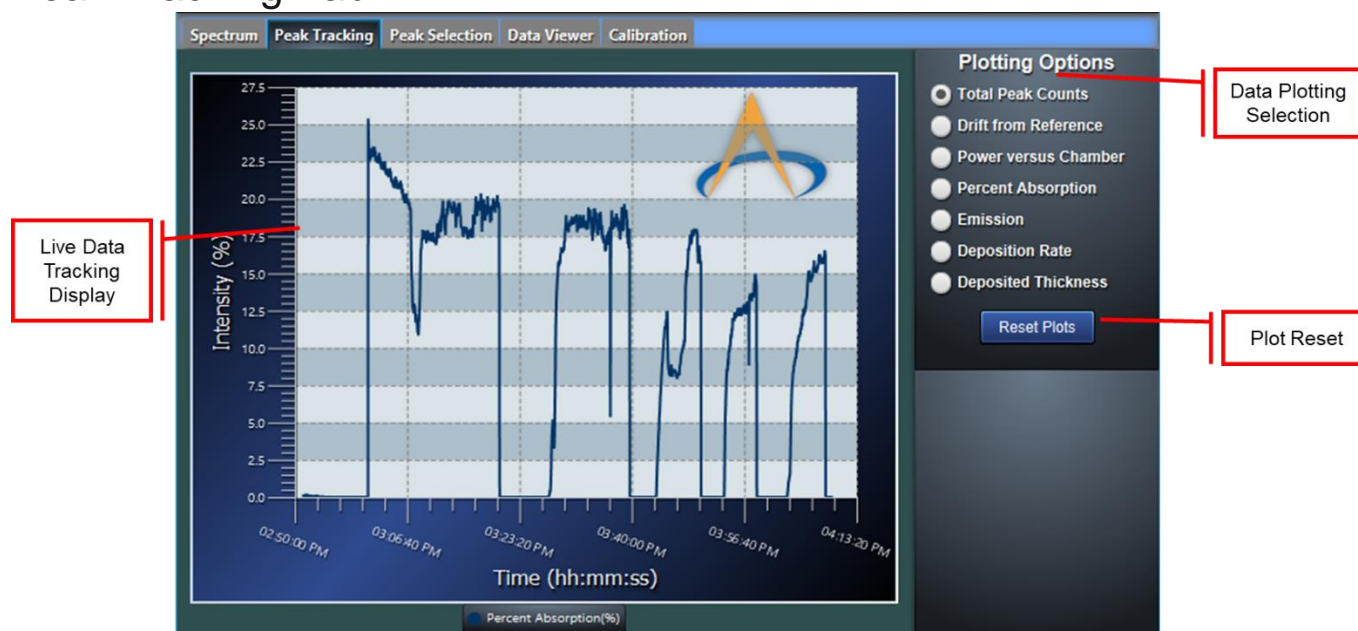

The peak tracking tab is the main tab for monitoring a process with the AtOMS system. The software provides a variety of ways to view the data. The panel includes data tracking options which can be selected from the list labeled **Plotting Options** on the right. The plotted data can be reset by using the “Plot Reset” button on the right side of the panel.

Drift from Reference

Tracks the difference between the total peak counts and the original reference peak.

Power versus Chamber

Tracks the difference between the elemental peak outside the chamber and inside the chamber.

Percent Absorption

Tracks the total percent absorption through the chamber. *This is the primary data that is used to track the deposition process.*

Emission

Tracks the total atomic emission occurring within the chamber during deposition.

Deposition Rate

Tracks the calculated deposition rate based on the calibration files provided for each element.

i Note

Requires calibration to display. Calibration is system dependent and individual instructions will be provided based on facility capabilities. Please contact AccuStrata for more information.




Deposition Thickness

Tracks the total cumulative thickness calculated from the beginning of the deposition.

i Note

Requires calibration to display. Calibration is system dependent and individual instructions will be provided based on facility capabilities. Please contact AccuStrata for more information.

Peak Selection Tab

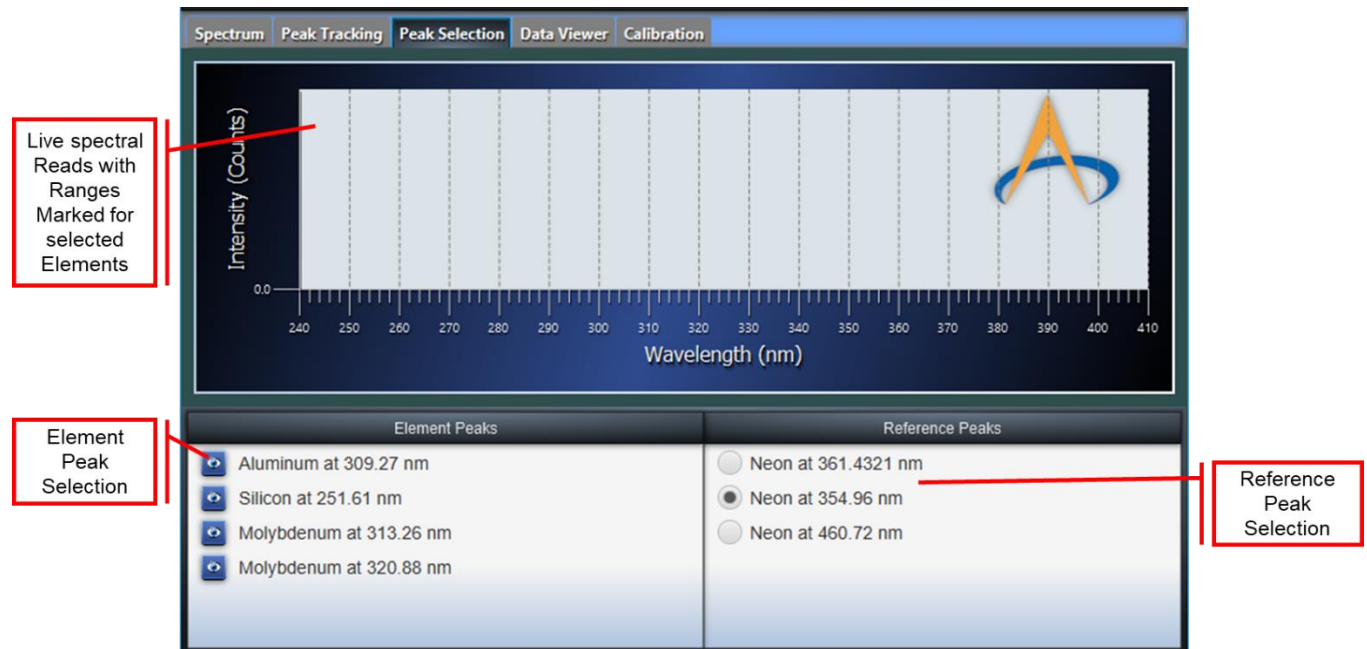

The Peak Selection Tab allows the user to select which elements to monitor as well as which reference peak to use. To change which elements are being monitored, toggle the “eyeball” next to each elemental line.


i Note

Settings on this panel can only be changed when a run is not occurring.

This panel will also display the live readings with range markers for each of the selected elemental peaks. The range indicated on the plot represents the region over which the peak is integrated.



Data Viewer Tab

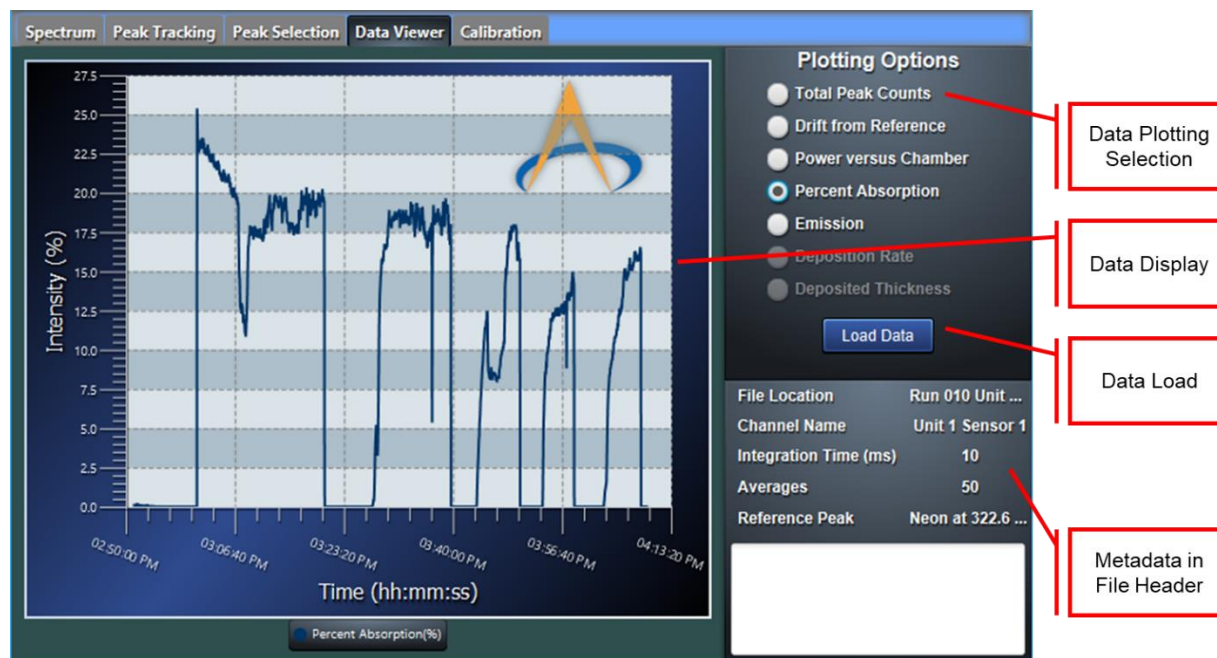


The Data Viewer Tab allows users to review old runs. The user can select between each of the plots by using the “Plotting Options” radio buttons. The metadata about the run including filename, sensor information, integration time, number of averages and the reference peak used for the data as well as any comments saved by the user will be displayed in the “Metadata” section. To load data, press the “Load Data” button and select the desired file.

WARNING

Only the processed files saved by the AtOMS system will display properly. Raw data files will not open successfully.



Calibration Tab

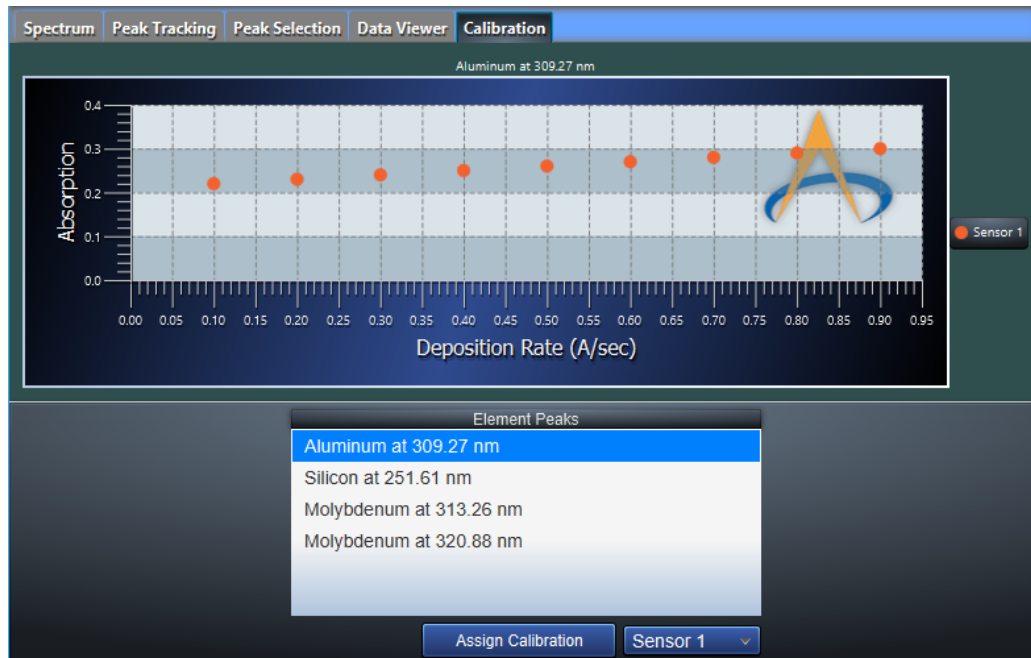

The Calibration tab is used to provide calibration curves that can be used to convert atomic absorption to deposition rate and thickness. To assign a calibration file to a specific elemental peak:

1. Selected the elemental peak that the file will be assigned to from the “Element Peaks” list.
2. Select the sensor that the calibration file if for from the dropdown menu.
3. Press the “Assign Calibration” button
4. Select the calibration file.

When an element is selected, the currently assigned calibration data for each sensor is displayed in the plot. If no calibration data is assigned, the plot will be blank.

To change a calibration file, carry out the assignment procedure and select the new file for the selected element.



System Menu

Setup Spectrometer Display

Set the display parameters for the spectrometer tab:

Sensors to Display

☒ Display Unit 1 Sensor 1

Spectra to Display

☐ Display Reference Spectrum ☐ Display Open Shutter Data ☐ Display Closed Shutter Data

Plot Limits

X Axis Limits

Minimum X: 300 Maximum X: 330

Y Axis Limits

Minimum Y: 0 Maximum Y: 10000

☐ AutoScale Data

OK Cancel

This menu allows changes to the Spectrometer tab plot.

Sensors to Display

Selects all sensors to be displayed on the Spectrum tab. Sometimes displaying all sensors at once can make viewing difficult. This option allows users to display only the spectra that are currently being monitor

Edit Peak Description

Peak Definition Utility

Available Peak Definitions

- ▶ Neon at 361.4321 nm
- ▶ Neon at 354.96 nm
- ▶ Neon at 460.72 nm
- ▶ Aluminum at 309.27 nm
- ▶ Silicon at 251.61 nm
- ▶ Molybdenum at 313.26 nm
- ▶ Molybdenum at 320.88 nm

Peak Definition Creator

Element: Aluminum

Central Wavelength (nm): 309.27

Minimum Wavelength (nm):

Maximum Wavelength (nm):

Is a Reference Peak: False

Buttons: Delete Select..., Add To List

Annotations:

- List of Currently Defined Peaks
- Add New Peak to List
- Delete Selection from List
- Element Selection
- Central Wavelength for the Element
- Lower Wavelength Value
- Upper Wavelength Value
- Define the Peak as a Reference or Elemental Peak

The Edit Peak Description menu allows users to create new peak definitions for various elements.

List of Currently Defined Peaks

A list of all currently defined peaks that will be selectable in the Peak Selection tab. Information about the element, wavelength range, central wavelength and whether the peak is considered a “references peak” can all be determined by expanding the tree list (down arrow).



Peak Definition Creator

To define a new peak to be used in the Peak Selection tab, fill in the values for the peak under this header. Every peak definition requires:

Element: The element that is being detected by the system. Only elements supported by AtOMS will be displayed in this list. If the desired element is not in the list, please contact customer support.

Central Wavelength: The elemental peak that will be monitored. Many elements have multiple elemental peaks that can be used to monitor. Most commercial HCL manufacturers will include the supported elemental peaks with each lamp.

Wavelength Range: Due to the precise nature of the atomic absorption process, many detectors are not perfectly calibrated for across the entire detector range. The “Minimum Wavelength(nm)” and “Maximum Wavelength” values are used to define the region in which the peak will be integrated. *These are user defined settings and can vary depending on element, detector and intensity.*


Many elements have multiple analytical lines. Most HCL manufacturers will include the most useful analytical lines on the HCL container. To the right is an example of an HCL label that comes with a Silicon HCL. The two recommended analytical lines are marked with red boxes. There may be other elemental lines not listed, only the most sensitive lines are included with the lamp container.



Note

Optimization of the minimum and maximum wavelength values is critical to system performance and should be tuned to provide maximum sensitivity and better signal to noise.

Hollow Cathode Settings

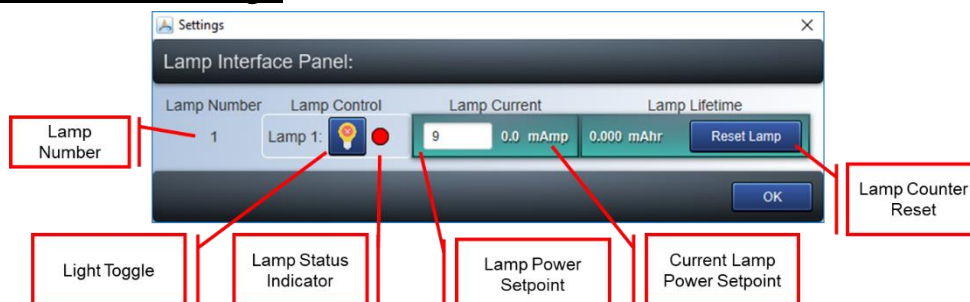



Lamp Toggled On

Indicates that the user intends for the lamp to be on.



Lamp Toggled Off

Indicates that the user intends for the lamp to be off.



System Status-Lamp On

The AtOMS system is providing power to the Hollow Cathode Lamp.




System Status-Lamp Off

The AtOMS system is not providing power to the Hollow Cathode Lamp.

Lamp Current

Provides users with the ability to set the HCL lamp current. Each HCL has different power settings.

To the right there is an example of a Silicon hollow cathode lamp package. Most lamps come with two power specifications. The first is the **Operating Current**. This is the current that is recommended by the manufacture for lamp operation. The second specification is the **Maximum Current**. This is the maximum current that the lamp can be set at without causing damage to the HCL. Both of these values are element and vendor specific, so it is critical to ensure that settings are entered correctly before operating the unit.

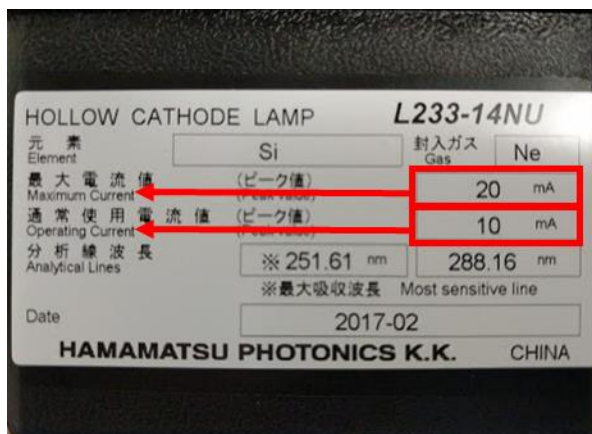


WARNING

Incorrect power settings for a lamp can lead to decreased lamp lifetime or catastrophic lamp failure. Use caution when setting up the lamp.




Detector Menu

Set Integration Time

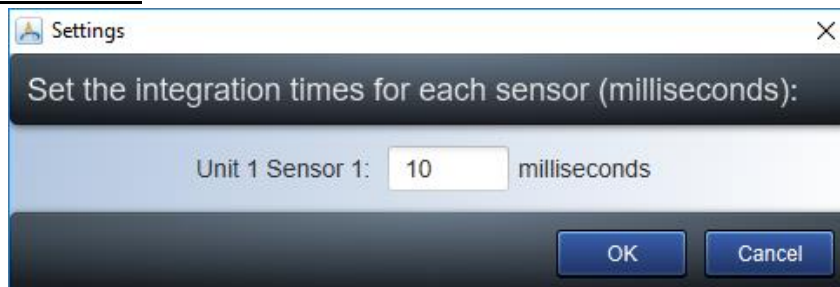


The **Integration Time** is the amount of time over which the detector will collect light. The higher the integration time, the higher the total peak counts will be. *The detector generally is most sensitive with counts around 40,000*; however, other factors such as reference peak intensity and lamp power may affect this.

Set Scans to Average

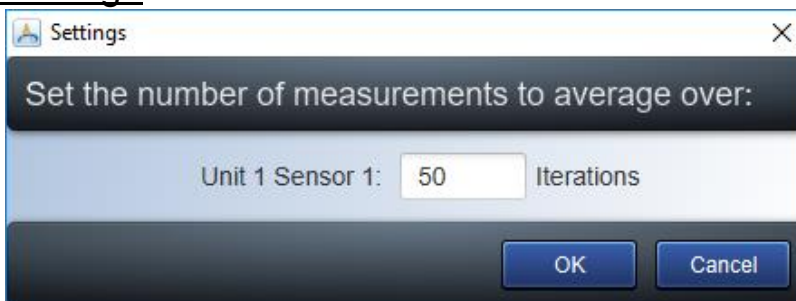

The **Scans to Average** is used to increase statistics for the measurement and improve the signal to noise; however, increasing the number of scans will also increase the data acquisition time. Users should balance the number of scans with the rate at which the processes are being monitored (sampling rate).

Set Shutter Ratio

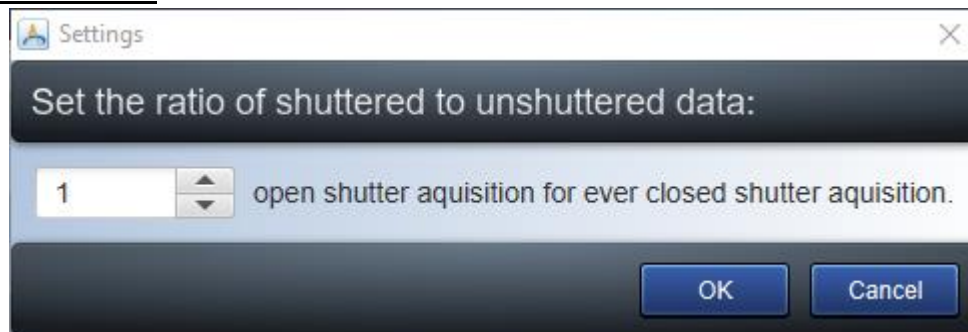

Shutter Ratio is the number of times a reference read will be reused. A setting of 1 will take a new reference reading for every measurement. This is the recommended setting and should only be changed by an advanced user.



Select Chambers



The AtOMS system can support up to 3 sensors. In some cases, not all sensors are needed. This menu option allows users to select which sensors will be measured. The channel number corresponds to the ports on the back of the AtOMS system.

Serial Communication

The AtOMS allows for communication via RS232 serial connector. The communication allows for control of the system as well as receiving calculated values from the system.


Hardware Setup


To configure the AtOMS hardware to support RS232 communicate, first plug the USB-to-RS232 converter into the AtOMS USB port on the front of the system. The system supports either USB port. The other USB port will be used for a keyboard/mouse pair. The RS232 can be removed at any time to accommodate the USB memory stick for file transfers.

Once the converter is plugged in, the correct COM port that was detected by windows must be determined. If it is unclear which



COM port is being used, right click on the Windows

button  on the bottom left and select "Device Manager". Under the list of hardware select "Ports (COM & LPT)". Here you will find the list of available COM port devices. There will be at least 2 COM devices. The "Arduino Mega 2560" is the AtOMS control board. The COM port labeled as USB-to-Serial Comm Port will be the COM port for the RS232.

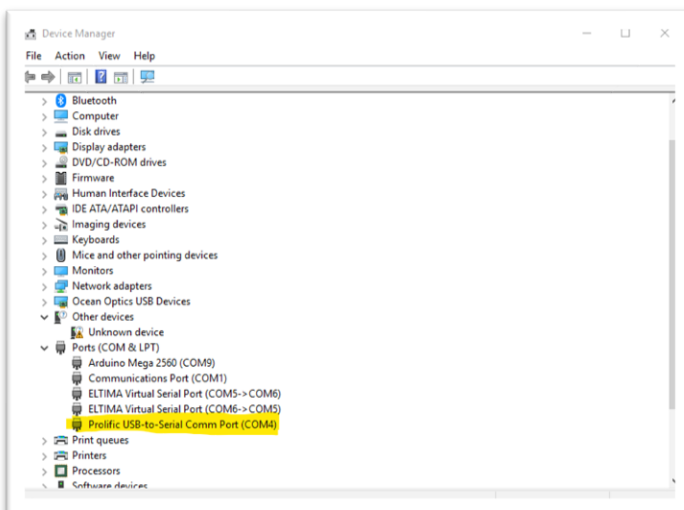




Software Setup

To configure the software for using the external communication, first go to “Communication Settings” under the “System” menu. This will bring up the settings panel. The settings panel will allow the user to set the Baud Rate and COM port for the external communication. The communication port list is auto-populated with the currently available ports, so the RS232 to USB converted must be plugged in prior to setting the COM port. The system assumes the following parameters:

Bits	8
Stop Bits	1
Parity	None

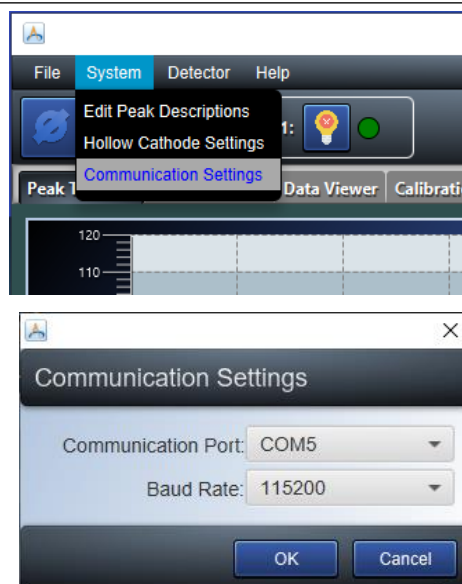

Communication Protocols

This section covers how to communicate with the AtOMS system and how to receive information from an external source. All communication is String based to provide a simple mechanism for communication.

Outputs

The system sends out a double value that is the last calculated absorption value followed by an carriage return and an end line. For example:

96.8/r/n

This would mean that the last Absorption reading was 96.8%.

Inputs

AtOMS provides external communication signals that can be used to control the status of the system. The following table lists all current communication capabilities:

Command	Parameter	Suffix	Effect	Button
Reference		/r/n	Take Reference	
Recording	(on/off)	/r/n	Toggle Recording	
Reading	(on/off)	/r/n	Toggle Reading	

For example:

Record on/r/n

Would start recording data. Stopping recording with “Recording off/r/n” will automatically save the data to the default location without any file suffix.



About



The About menu is used to identify critical versioning information about the software. To ensure that updates have been correctly implemented, use the Version number.

Contact

Support:

AccuStrata Incorporated
11900 Parklawn Dr., Suite 203
(240) 223-5400
support@accustrata.com